

SLH • SLS • SLHS

Лабораторные плитки/мешалки

SI Analytics



По вопросам продаж и поддержки обращайтесь:

Архангельск (8182)63-90-72
Астана +7(7172)727-132
Белгород (4722)40-23-64
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89
Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Казань (843)206-01-48

Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Липецк (4742)52-20-81
Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41

Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Новосибирск (383)227-86-73
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78

Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Тверь (4822)63-31-35
Томск (3822)98-41-53
Тула (4872)74-02-29
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Ярославль (4852)69-52-93

сайт: <http://sianalytics.nt-rt.ru> || эл. почта: ssc@nt-rt.ru

SLH, SLS, SLHS – new laboratory heatplates and stirrers from SI Analytics

New laboratory equipment from SI Analytics: the SLH laboratory heatplate, SLS stirrer and the SLHS heatplate/stirrer combination.

The **compact design** with a footprint of 205 x 260 mm allows the use on crowded laboratory tables or under fume hoods. Yet, the working surface of 190 x 190 mm makes the SLH, SLS and SLHS compatible with most standard size borosilicate bottles, beakers and flasks.

All three units feature a **ceramic coated stainless steel plate** highly resistant against strong acids and bases. The powder coated aluminium cast housing is heat resistant and can withstand heavy laboratory use.

The stirrer is equipped with a **speed control knob for a range of 60–500 rpm**. The drive motor is equipped with ball bearings for smooth running and a long service life.

The heatplate of the SLH and SLHS has **maximum power consumption of 500 Watts** and is electronically controlled to prevent overcharging.

Two control LEDs on the front plate light up when heating and stirring function are on.

Technical Data

Model	SLS-110 and SLS-230	SLH-110 and SLS-230	SLHS-110 and SLHS-230
Type	Laboratory stirrer	Laboratory heatplate	Laboratory stirrer/heatplate
Stirrer speed range	60–1500 U/min	N/A	60–1500 U/min
Temperature range	N/A	Room temperature +5°C–380 °C	
Control unit	Breakerless electronic control		
Material (stirring/heating plate)	Ceramic coated stainless steel		
Physical dimensions	Plate	190 x 190 mm / 7,5 x 7,5 “	
	Instrument	205 mm width x 260 mm depth x 110 mm height / 8,5 x 11 x 4,5”	
Power input	Max. 15 Watts, 1.0 A	Max. 500 Watts, 3.0 A	
Voltage	110V, 60Hz or 230V, 50/60 Hz		

Ordering information

Product	Description	Type No.	Order No.
Laborheizrührer SLHS	Ceramic coated stainless steel plate, 190x190 mm, 60–1500 rpm, 0,5 kW heating power, with mains cable, 230 V	SLHS-230	285416400
Laborheizrührer SLHS	Ceramic coated stainless steel plate, 190x190 mm, 60–1500 rpm, 0,5 kW heating power, with mains cable, 110 V	SLHS-110	285416410
Laborheizplatte SLH	Ceramic coated stainless steel plate, 190x190 mm, 0,5 kW heating power, with mains cable, 230 V	SLH-230	285416420
Laborheizplatte SLH	Ceramic coated stainless steel plate, 190x190 mm, 0,5 kW heating power, with mains cable, 110 V	SLH-110	285416430
Laborrührer SLS	Ceramic coated stainless steel plate, 190x190 mm, 60–1500 rpm, with mains cable, 230 V	SLS-230	285416440
Laborrührer SLS	Ceramic coated stainless steel plate, 190x190 mm, 60–1500 rpm, with mains cable, 110 V	SLS-110	285416450

По вопросам продаж и поддержки обращайтесь:

Архангельск (8182)63-90-72
Астана +7(7172)727-132
Белгород (4722)40-23-64
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89
Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Казань (843)206-01-48

Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Липецк (4742)52-20-81
Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41

Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Новосибирск (383)227-86-73
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78

Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Тверь (4822)63-31-35
Томск (3822)98-41-53
Тула (4872)74-02-29
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Ярославль (4852)69-52-93